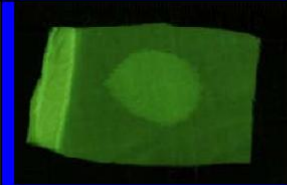

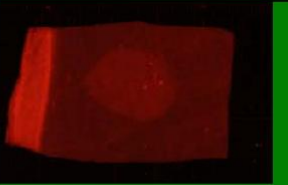
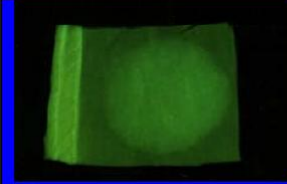
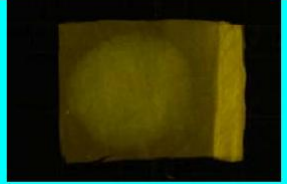

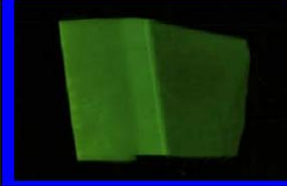

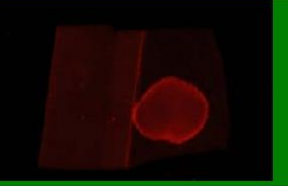


Locating Flammable Liquids Using the Crime-lite Range

It is common to find traces of accelerant at arson scenes and in other insurance cases investigations where a fire has been involved. Identification of an accelerant provides information about where and how the fire started. It may also give an indication whether the fire was accidental or started with intent.

Flammable liquids were applied to a white cloth and allowed to dry. The flammable liquids then gave fluorescent emissions when illuminated using the Crime-Lite range and viewed and photographed with their respective filters.

The images shown were photographed using the DCS4 system and the Crime-Lite 8x4 using the new anti-glare viewing filters.

Accelerant	Blue (445nm) With GG495 Filter	Blue/Green (480nm) With OG550 Filter	Green (530nm) With OG590 Filter
Petrol			
Diesel			
Methylated Spirits			

Sales Support and Feedback - Foster + Freeman welcome feedback from Customers regarding this product. Please contact one of our offices if you would like to pass on your comments. Foster + Freeman are pleased to offer advice, installation, training and on-site maintenance worldwide for all of their products.

FOSTER + FREEMAN LTD
 25 Swan Lane
 Evesham
 Worcestershire
 WR11 4PE
 UK
 ☎ + 44 (0) 1386 41061
 ☎ + 44 (0) 1386 765351
 sales@fosterfreeman.com
 www.fosterfreeman.com

Head Office & UK Sales Office

Foster + Freeman reserve the right to alter the specification of this product, accessories and consumables without prior notice.
 This document contains proprietary information that is protected by copyright.
 All rights are reserved.
 No part of this publication may be reproduced in any form whatsoever without the prior, written permission of Foster + Freeman Ltd.
 Copyright © Foster + Freeman Ltd

FOSTER + FREEMAN
 USA INC
 46030 Manekin Plaza
 Sterling
 Virginia 20166
 USA
 ☎ +1 888 445 5048
 ☎ +1 888 445 5049
 usoffice@fosterfreeman.com
 www.fosterfreeman.com

US Sales Office